



*Your best portable
sample preparation
partner*

Tissue Grinders

Features

- Handheld, motor-driven
 - Re-suspends pellets or disrupts soft tissue in microcentrifuge tubes
 - Easy to operate
- Avoid Contamination
 - Disposable plastic grinding pestle
- Battery Powered
 - Can be used in mobile applications
 - Ideal for field testing
- Easy to clean
 - Stainless and plastic grinding pestles can be autoclaved
- Multiple Applications
 - Plant and animal tissues; DNA/RNA extraction

Molecular Diagnostics For Everyone

Labguide Co., Ltd. (Taiwan)

E-mail : terry@seed.net.tw

Tel. : +886-2-29406162

Add. : 3F., No.37, Ln. 232, Huaxin St., Zhonghe Dist.,
New Taipei City 235, Taiwan (R.O.C.)

<http://www.labguide.com.tw>

Labguide Co., Ltd. (Shenzhen)

E-mail : admin@labguide.com.cn

Tel. : +86-075527750787

Add. : Address: C603, Baohe Building, No.98 of Baonong
North Alley, Bao'an District, Shenzhen, China.

<http://www.labguide.com.cn>













Animal Tissue Samples

Samples	krpm	Processing Time
Heart	10	High
Liver	5	Low
Spleen	5	Medium
Lung	8	Medium
Kidney	8	Medium
Small intestine	9 - 10	High
Large intestine	5	Low
Brain	5	Low
Testis	8	Medium
Ear	10	N/A
Skin	10	N/A
Marrow	5	Low
Tonsil	8	Medium
Thyroid	8	Medium
Lymph	10	High
Dorsal muscle	10	Medium
Vessel tissue	10	N/A


Plant Tissue Samples

Samples	krpm	Processing Time
Soybean leaf	6	Low
Wheat leaf	6	Low
Tobacco leaf	6	Low
Chinese Wildrye leaf	5	Low
Bamboo leaf	6	Low
Rice leaf	6	Low
Rice root	7	High
Rice stem	10	N/A
Rice seed	10	N/A
Arabidopsis thaliana seed	7	High
Arabidopsis thaliana leaf	6	Low
Arabidopsis thaliana root	5	Low

Performance

	Before	After
Tobacco leaf		
Lung		
Large intestine		
Wheat leaf		
Liver		
Heart		

Specifications

Model:	G50	G10
Motor		Large torque brushless motor with 50W constant power output
Speed	Adjustable speed ranging from 3000 - 8000 rpm. Speed shown on LCD screen	Large torque DC motor, low noise, with constant speed
Grinding pestle (Stainless/Plastic)	Grinding pestles can be autoclaved	12000 rpm
Power Supply	12V DC universal power supply, 12V DC battery pack or car charged	Grinding pestles can be autoclaved
Dimensions	170*50*26 mm	3.7V rechargeable lithium battery, up to 10 hrs of operation
		155*φ25 mm

Covate Bioscience Co., Ltd.

Covate Bioscience USA Inc.

Labguide Co., Ltd. (Taiwan)

E-mail : terry@seed.net.tw

Tel. : +886-2-29406162

Add. : 3F., No.37, Ln. 232, Huaxin St., Zhonghe Dist., New Taipei City 235, Taiwan (R.O.C.)

<http://www.labguide.com.tw>

Labguide Co., Ltd. (Shenzhen)

E-mail : admin@labguide.com.cn

Tel. : +86-075527750787

Add. : Address: C603, Baohe Building, No.98 of Baonong North Alley, Bao'an District, Shenzhen, China.

<http://www.labguide.com.cn>